

INSTALLATION INSTRUCTIONS - PLUMBING

Fume Cupboard Plumbing Connections:

1. All the services are available on the side of the fume cupboard (max. 6 on each side).
2. Mains access needs to be max. 100mm from side wall position. The services side left/right was defined during quoting process and needs to be followed.
3. S/S braided hoses 1100mm long fitted with 1/4" BSP female connectors permit the connection to the mains. Each hose is labelled identifying its intended use.

School Fume Cupboard Plumbing Connection:

1. All the services are available on the bottom bench of the fume cupboard (max. 4 mounted on the bench)
2. S/S braided hoses 1100mm long fitted with 1/2" BSP female connectors permit the connection to the mains. Each hose is labelled identifying its intended use.
3. Connection:
 - a. Local delivery – fitted by Dynaflow.
 - The controls will be installed to the bottom bench by Dynaflow installation personnel.
 - The site plumber needs to connect the line to the mains using the supplied hose.
 - b. Remote installations – fitted by site plumber.
 - The controls will be installed to the bottom bench by the site plumber. Parts are shipped separately to avoid scratches.
 - The site plumber connects the service to the mains using the supplied components all the fittings are 1/2" BSP



Accredited No. 2910

Quality
Endorsed
Company

Errington Site
ISO 9001 LIC QEC23606 SAI Global Dynaflow Pty Ltd A.B.N. 28 008 493 178

30 Waratah Street
Errington NSW 2115
Australia

Phone: +61 (0) 2 9858 0190
Fax: +61 (0) 2 9858 1177

153 Magill Road
Stepney SA 5069
Australia

Phone: +61 (0) 8 8362 9099
Fax: +61 (0) 8 8362 9622